


Mounting H122SZX Focus Motor on Olympus SZX



1. Remove screw holding focus knob assembly.

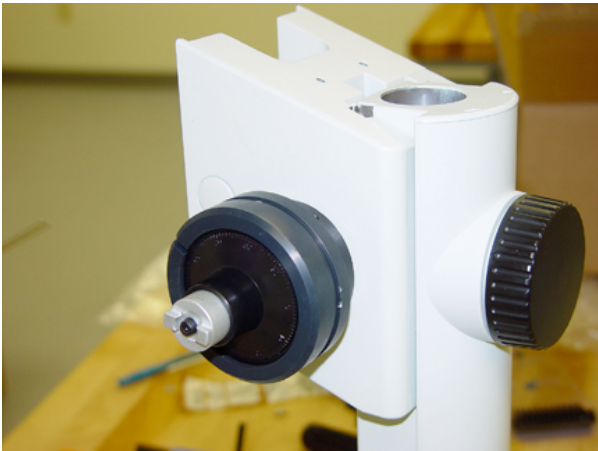

2. Attach hub to assembly using M3 x 25 socket screw and washer provided with focus motor assembly. Take care to make sure the hub and fine focus knob are concentric.



3. Slide the split ring sleeve over the coarse knob. The chamfered end of the sleeve should be outward from the body of the microscope.



4. Slide the focus motor bell housing over the sleeve. Tighten the 3 set screws on the outer diameter of the bell housing. This will clamp the housing to the coarse knob outer diameter



5. Mount the 5mm hub and black anti backlash coupler to the end of the focus motor. The hub will tighten to the motor shaft via the set screw on its side.



6. Slide the focus motor into the bell housing. While applying gentle pressure inward on the motor, tighten the thumb screw. Note you may have to rotate the motor in order to seat the coupling correctly.

End of Instructions

17 September 2008