


INSTRUCTIONS FOR MOUNTING Z ENCODER PROBE BRACKET
MODEL H394 TO LEICA DMIR INVERTED MICROSCOPE


1. Mount encoder reference bracket onto nosepiece as shown below. Note you will need to remove the glass contact plate from the rounded end of the assembly.

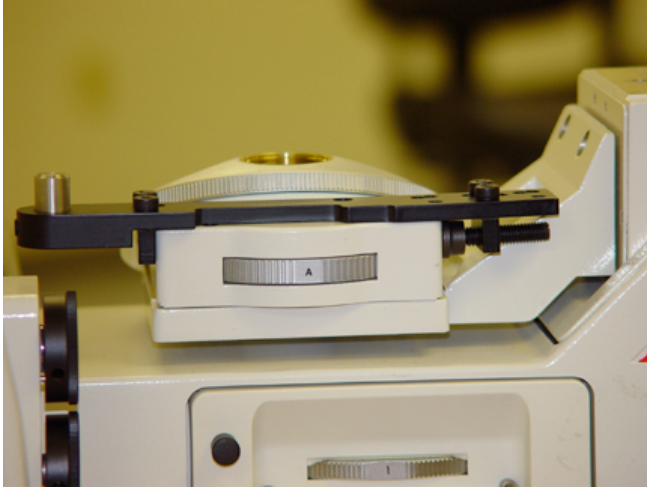

2. Mount the rod and probe clamp assembly onto the probe reference bracket as shown below. The rod and probe clamp assembly may be found loose in your kit of parts or it may already be installed onto the probe main bar.



3. Install Encoder probe into probe clamp and tighten clamp as shown below.

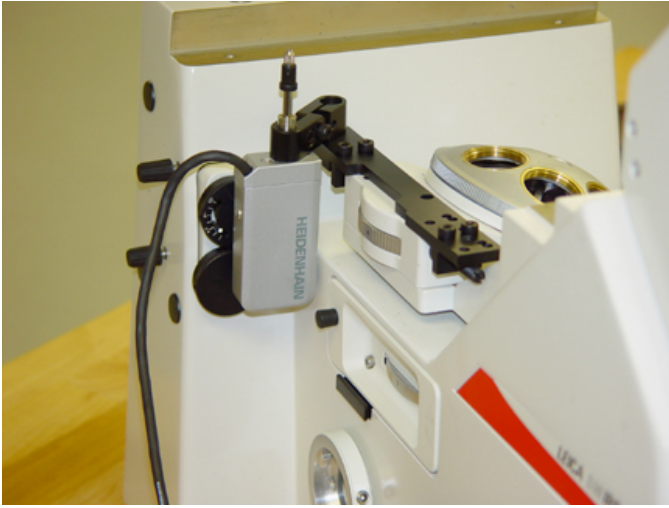

4. Mount stage onto microscope. Ensure that probe maintains contact with stage bottom throughout the required range of motion. The height of the encoder can be adjusted by moving it up or down within the probe clamp.



End of Instructions – JH 6/26/03