



## Z Encoder Probe Bracket Mounting Instructions – TE2000


1. Remove (2) M5 Nosepiece mounting screws

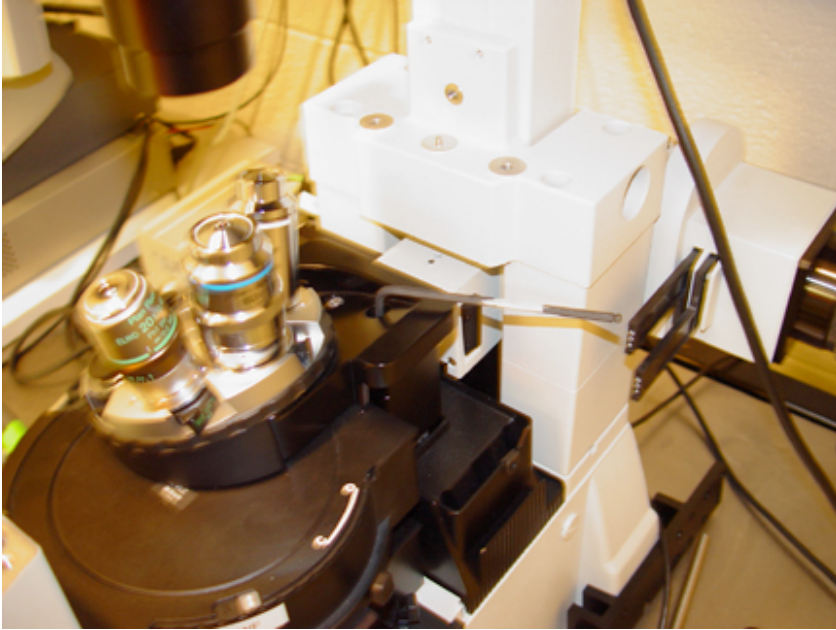


2. Place probe bracket on top of nosepiece and screw bracket thru the top of nosepiece and into the microscope using the 2 M5 x 40 captive fasteners supplied.

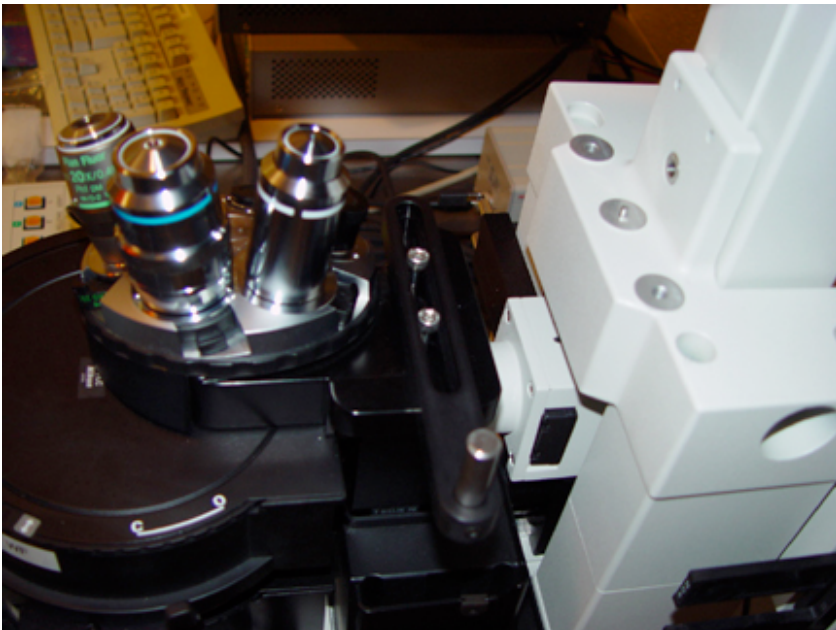

3. Mount the encoder probe to the stainless rod using the mounting clamps provided.

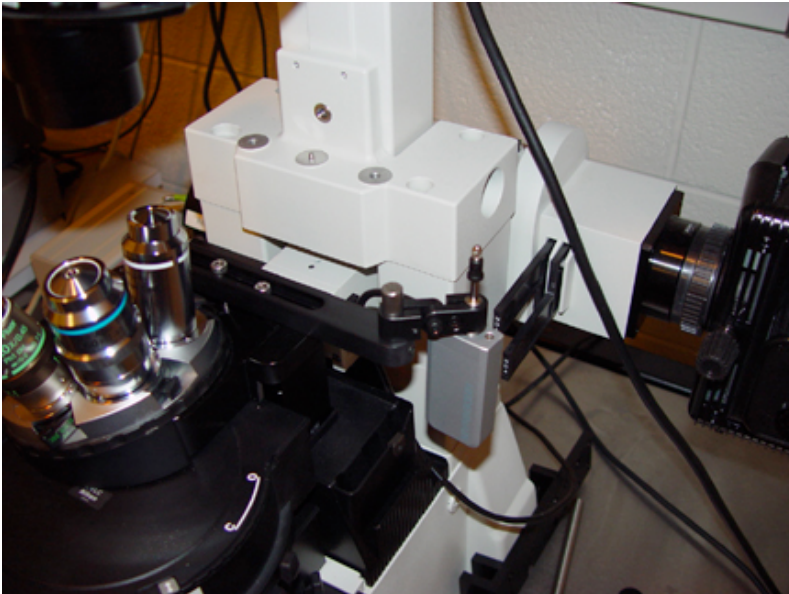

4. Install the stage and adjust the position of the encoder probe. Adjust the XY position of the bracket using the rotation of the mounting clamps and the slot in the mounting bracket so that the probe contacts the flat part of the bottom plate of the stage. Adjust the probe up and down on the stainless rod so that the encoder maintains contact with the stage bottom plate over the required Z travel.

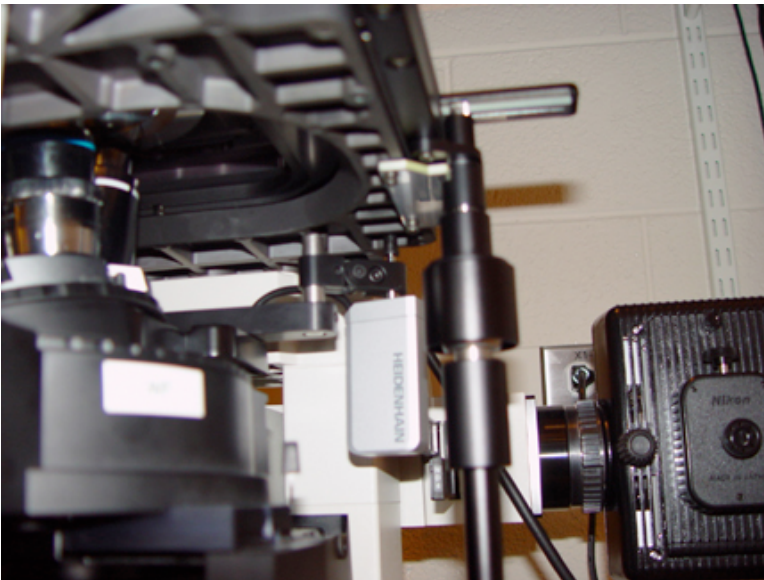

End of Instructions – JH 6/28/04