



IMPORTANT SAFETY INFORMATION



OPTISCAN III



PLEASE READ THE MANUAL BEFORE USING!



THE MANUAL IS AVAILABLE FROM PRIOR SCIENTIFIC AND SHOULD BE READ AND UNDERSTOOD BEFORE INSTALLATION AND USE. THE MANUAL IS AVAILABLE FROM WWW.PRIOR.COM

- Adhere to all warnings on the product and in the user manual.
- This product **MUST NOT** be exposed to open flames, extremely hot or cold temperature, water or moisture. **DO NOT** allow objects to fall, or liquids to spill, on the product.
- The OptiScan III is a Class I system and must be earthed.
- Connect the AC power cord only to designated power sources as marked on the product. Make sure the electrical cord is located so that it will not be subject to damage.
- Ensure the connection to the power source and the front panel power switch are easily accessible.
- Do not connect or disconnect anything to or from the ProScan III unless the controller is turned off first.
- **DANGER – Never alter the AC cord or plug.** The power cord set must be an appropriately rated and approved cord set in accordance in the regulations of the country it is used in. If the supplied plug adapter is not the correct fitting for your geographic area or if you are unsure about the relevant regulations, please contact your supplier for advice.
- **Do not attempt to disassemble, tamper or repair this product; it contains NO user serviceable parts and doing so may void the warranty.** Service should be performed **ONLY** by authorised service centres.
- Clean only the exterior of this product using a damp, lint free cloth.
- **Please contact Prior or an approved representative if you have any questions or comments.**



ALWAYS MAKE SURE THE POWER TO THE OPTISCAN III IS OFF BEFORE CONNECTING OR DISCONNECTING ANY COMPONENTS




OPTISCAN III JOYSTICK

Forward/backward - Y axis movement
Left/Right - X axis movement

Press to alter speed of focus motor

Press to alter speed of stage

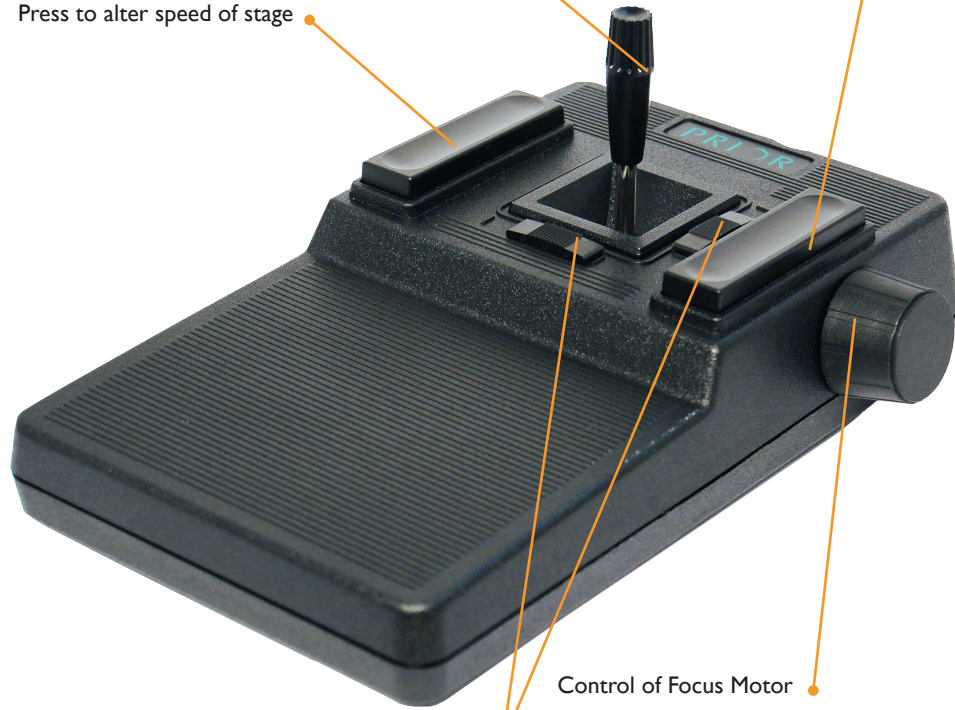

Control of Focus Motor

Tensioners - latch to hold joystick in off position

OPTISCAN III CONNECTIONS

RS232-1 Connection

RS232-2 Connection



Power

Joystick

USB

Z- Axis

XY Stage

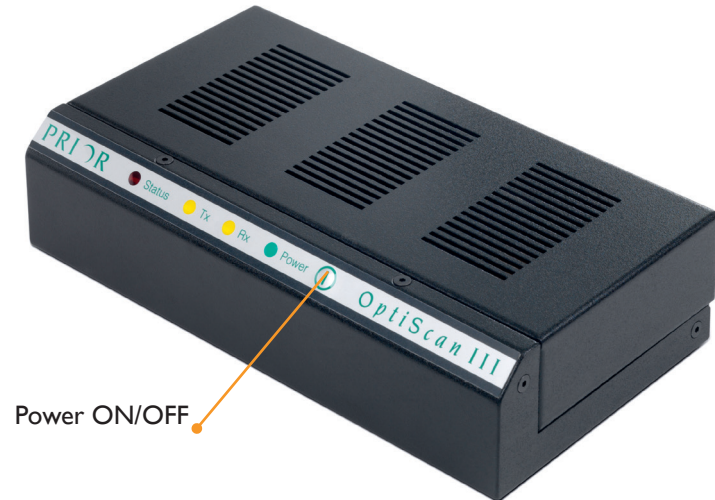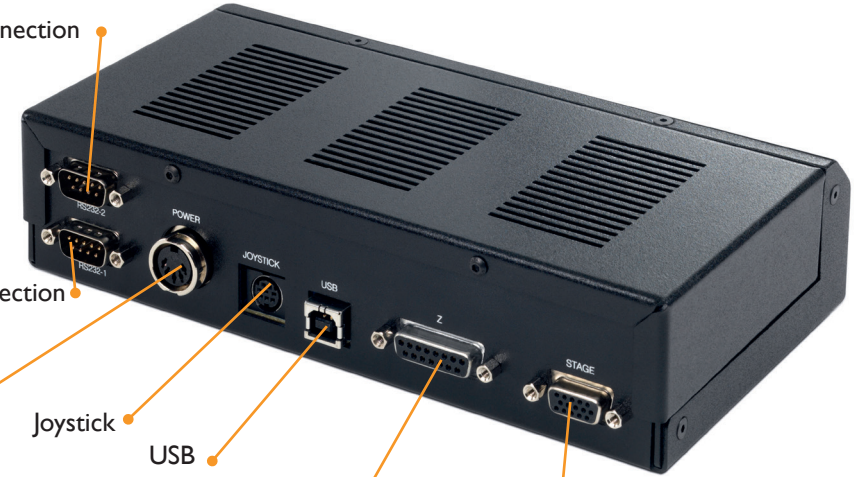

ALWAYS MAKE SURE THE POWER TO THE OPTISCAN III IS OFF BEFORE CONNECTING OR DISCONNECTING ANY COMPONENTS!

READ THE MANUAL BEFORE USE (WWW.PRIOR.COM)

