



## IMPORTANT SAFETY INFORMATION



### PLEASE READ THE MANUAL BEFORE USING!

THE MANUAL IS AVAILABLE FROM PRIOR SCIENTIFIC AND SHOULD BE READ AND UNDERSTOOD BEFORE INSTALLATION AND USE.


THE MANUAL IS AVAILABLE AT [WWW.PRIOR.COM/MANUALS](http://WWW.PRIOR.COM/MANUALS)

THE SOFTWARE IS AVAILABLE AT [WWW.PRIOR.COM/SOFTWARE](http://WWW.PRIOR.COM/SOFTWARE)

- Adhere to all warnings on the product and in the user manual.
- This product MUST NOT be exposed to open flames, extremely hot or cold temperature, water or moisture. DO NOT allow objects to fall, or liquids to spill, on the product.
- The ProScan III is a Class 1 System and must be earthed.
- Connect the AC power cord only to designated power sources as marked on the product. Make sure the electrical cord is located so that it will not be subject to damage.
- Ensure the connection to the power source and the front panel power switch are easily accessible.
- Do not connect or disconnect anything to or from the ProScan III unless the controller is turned off first.
- **DANGER – Never alter the AC cord or plug.** The power cord set must be an appropriately rated and approved cord set in accordance in the regulations of the country it is used in. If the supplied plug adapter is not the correct fitting for your geographic area or if you are unsure about the relevant regulations, please contact your supplier for advice.
- Do not attempt to disassemble, tamper or repair this product; it contains no user serviceable parts and doing so may void the warranty. Service should be performed ONLY by authorised service centres.
- Clean only the exterior of this product using a damp, lint free cloth.
- Please contact Prior or an approved representative if you have any questions or comments.



## ProScan III Connections

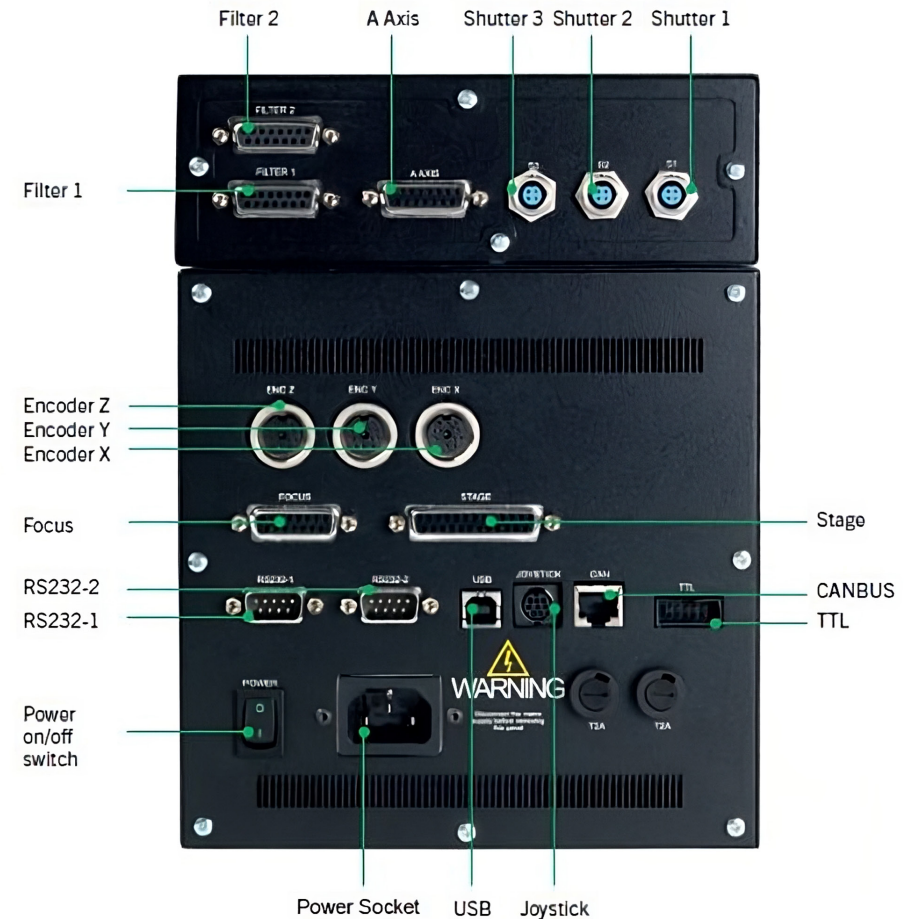


The ProScan III utilizes automatically installing Windows-compatible drivers.

**DO NOT CONNECT OR DISCONNECT ANYTHING TO OR FROM THE PROSCAN III UNLESS THE CONTROLLER IS TURNED OFF FIRST.**




# ProScan III Connections


## Control

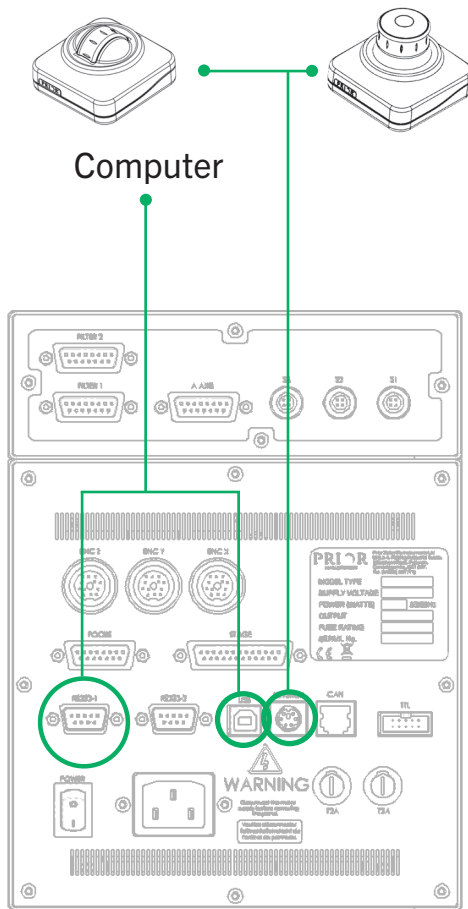


## Stage

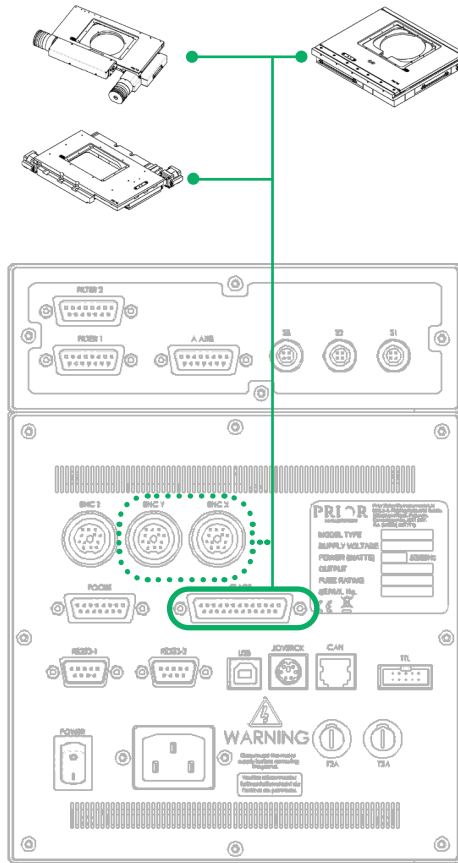

## Focus

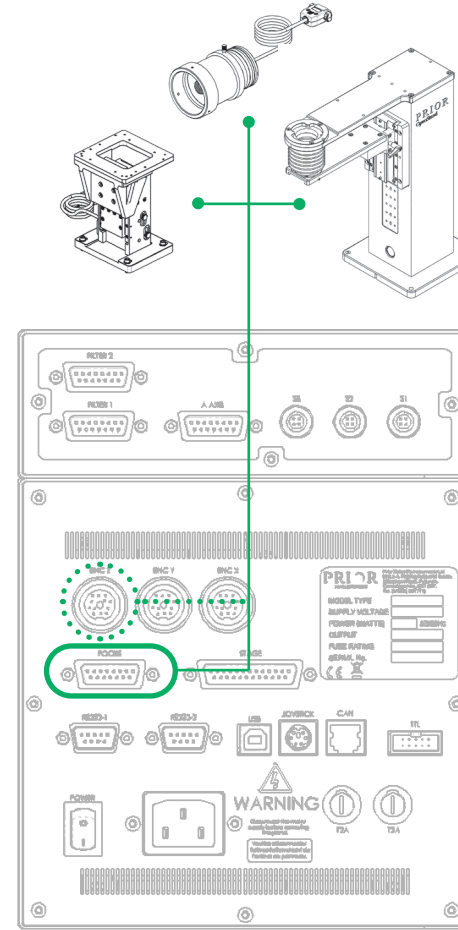

## Illumination

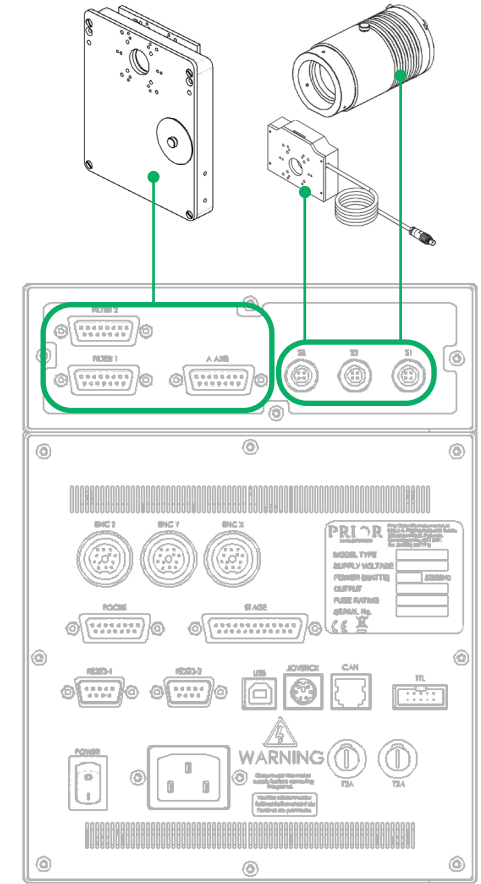

- Connection between ProScan III and Prior Scientific device
- ..... Encoded connection between ProScan III and encoded Prior Scientific device

- Connection between ProScan III and Prior Scientific device
- ..... Encoded connection between ProScan III and encoded Prior Scientific device

**DO NOT CONNECT OR DISCONNECT ANYTHING TO OR FROM THE PROSCAN III UNLESS THE CONTROLLER IS TURNED OFF FIRST.**

**ENSURE THE CORRECT ENCODER CABLE GOES INTO THE CORRECT PORT READ THE MANUAL BEFORE USE (WWW.PRIOR.COM/MANUALS)**



ProScan III Quick Start Guide © Prior Scientific 2023  
W4641 V4



Visit [www.prior.com](http://www.prior.com) for more information & download the software from [www.prior.com/software](http://www.prior.com/software)