

How to reset a Prior Controller

RESETTING YOUR CONTROLLER

The ES10ZE, OptiScan and ProScan range of controllers can be reset to their default settings with these instructions. **Note that resetting the controller will remove all settings from the controller and return them to default, so ensure you note down any settings you wish to keep.**

The reset command is sent via PRIOR TERMINAL software. For help with installation of Prior Terminal see the guide 'Installing and Using Prior Software'.

The command can be sent via RS232 port (if there is more than one on the controller, it doesn't matter which one is connected to the computer) or USB port.

In all cases you should send the command RESET. Press the 'Enter' key to enter the command. The controller will respond with '0' and possibly with 'CODE UPDATED – INITIALISE'. This will signify your controller has been reset and is now ready to use.

If using a USB connection the connection may cease. This no cause for alarm; simply restart the controller.